

BSC Failure or Alarm

Remain Calm



Immediately Stop Work



Close All Primary Containers

if not working with biohazardous materials, surface decontaminate items, remove them from BSC, post “Do Not Use” signage and report problem to supervisor; if working with biohazardous materials continue to follow the instructions below



Remove Gloves and Discard in BSC Waste



Turn Off BSC and Close Sash



Wash Hands



Post “Do Not Use” Signage



After 30 Minutes, Don Clean PPE, Surface Decontaminate Then Remove Items in BSC



Decontaminate Work Surface, Close Sash, Leave “Do Not Use” Signage on BSC



Report Problem to Supervisor